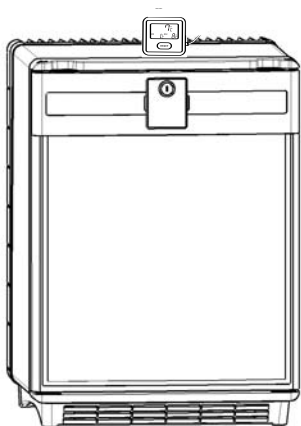
Dometic miniCool

DS301H DS601H
Type: DSH20-60

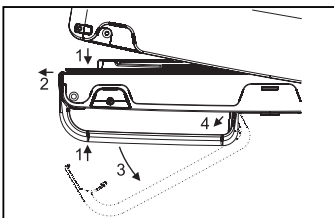
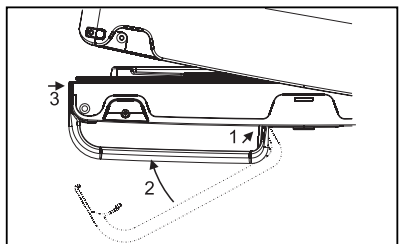
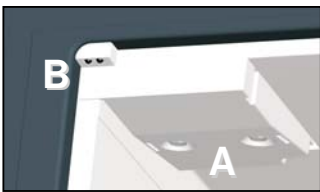
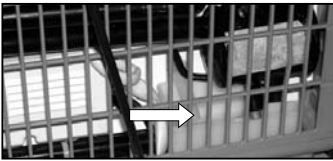
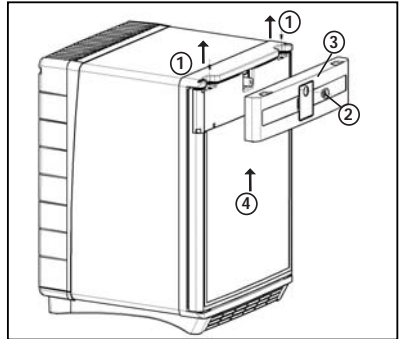
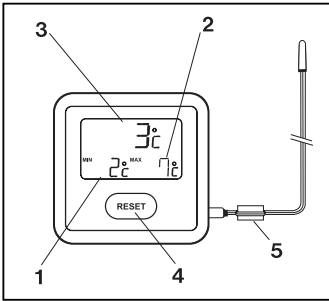
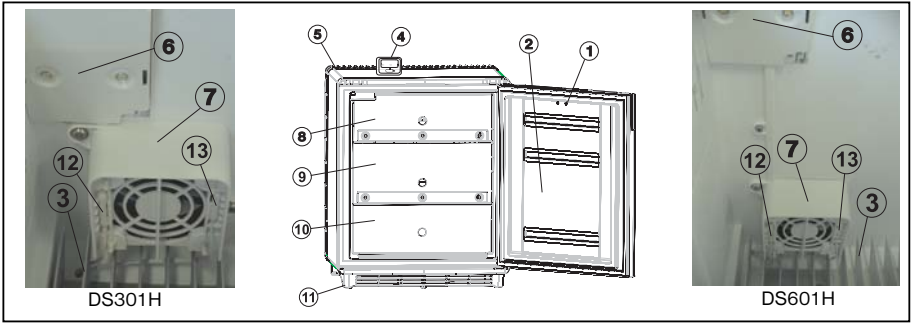
Health Care refrigerator
Arzneimittel-Kühlgerät
Réfrigérateur pour Soins de santé
Frigorifero medicale
Koelkast voor medische doeleinden
Refrigeração medicinal
Kylskåp för läkemedel
Medisinkjøleskap
Køleboks til sundhedssektoren
Näyte/lääkejääkaappi
Frigorífico Health Care
Kórházi hűtőszekrény

Operating Instructions
Bedienungsanleitung
Mode d'emploi
Istruzioni per l'uso
Handleiding
Käyttöohje
Bruksanvisning
Bruksanvisning
Betjeningsvejledning
Manual de instrucciones
Kezelési tájékoztató

GB
DE
BE
IT
NL
SF
SE
NO
DK
E
HU

Please read these operating instructions carefully before putting the refrigeration unit into operation. If you later sell or dispose of it, please ensure that the new owner receives these operating instructions.

Please follow the installation details below carefully. Guarantee is valid for products installed as described only.

Thank you for choosing our appliance. We are sure it will provide you with trouble-free use.

In the following, we would like to familiarise you with some symbols, which we bring to your attention to ensure the safe and efficient operation of the appliance:

source of danger, in event of improper operation

suggested useful tips to read

information concerning environmental protection

This refrigerator is intended for the storage of drugs at temperatures within the range +2°C to +8°C. A digital type thermometer is provided which may be fixed to the top of refrigerator. (Fig. 1., page 4)

Attention!

THIS REFRIGERATOR IS NOT SUITABLE FOR THE STORAGE OF BLOOD.

Attention!

• **This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience and knowledge, unless they have been given supervision or instruction concerning use of the appliance by a person responsible for their safety.**

Children should be supervised to ensure that they do not play with the appliance.

- Any electrical work required to install this appliance should be carried out by a qualified electrician.
- This product should be serviced by an authorised Service Engineer, and only genuine spare parts should be used.
- It is dangerous to alter the specifications or modify this product in any way.
- Care must be taken to ensure that the appliance does not stand on the electrical supply cable.
- There are working parts in this product which heat up. Always ensure that there is adequate ventilation as a failure to do this will result in component failure and possible content loss. See installation instructions.
- Parts which heat up should not be exposed. The back of the product should be close to a wall but leaving the required distance for ventilation as stated in the installation instructions.
- Before defrosting, cleaning or maintenance work is carried out, be sure to switch off the appliance and unplug it.
- **DO NOT USE SHARP INSTRUMENTS** to scrape off frost or ice. Under no circumstances should solid ice be forced off the cooling fins. Solid ice should be allowed to thaw when defrosting the appliance. See defrost instructions.
- This appliance is heavy. Care should be taken when moving it.
- Under no circumstances should you attempt to repair the appliance yourself. Repairs carried out by inexperienced persons may cause injury or more serious malfunctioning. Refer to your local Service Centre and always insist on genuine spare parts.
- In this appliance the storage of any toxic or explosive substance is forbidden!

CONTENTS

1. Unpacking
2. View of the appliance
3. Cleaning
4. Positioning the appliance
 - 4.1. Installation instruction
5. Using the appliance
 - 5.1. Initial operation
 - 5.2. Temperature regulation
6. Thermometer
7. Useful suggestions
8. Defrosting, cleaning and maintenance
9. Interior light
10. Replacing the decorative door panel
11. Fitting up and removing the handle (optional)
 - 11.1. Fitting up the handle
 - 11.2. Removing the handle
12. Customer service
13. Putting into operation, technical data
 - 13.1. Connecting to electrical mains
14. Environmental protection information
15. Recycling

**Attention!**

Warranty arrangements are in accordance with EC Directive 44/1999/CE and the normal conditions applicable in the country concerned.

For warranty or other servicing, such as spare-parts, please contact our Dometic Service Network.

The warranty does not cover any damage due to improper use.

The warranty does not cover any modifications to the appliance or the use of non-original Dometic spare-parts.

The warranty does not apply if the installing and operating instructions are not adhered to.

When contacting Dometic Service Network, please state the model, product number and serial number. You will find this information on the data plate inside of the refrigerator.

1. Unpacking

After removal from the cardboard packaging, make sure the appliance is not damaged. If you find damage to the

appliance resulting from transport, report it immediately to the transportation firm.

2. View of the appliance (Fig 1)

- | | |
|----------------------------|--|
| 1. Door lock | 8. Upper tray (Cutout on Upper Left Corner of Front Plate) |
| 2. Door | 9. Middle tray (With Highest Front Plate) |
| 3. Evaporator | 10. Lower tray (Injection-moulded Compact Component) |
| 4. Thermometer | 11. Footframe |
| 5. Cooling unit room cover | 12. Sensor of Fuzzy Logic |
| 6. Interior light | 13. Sensor of thermometer |
| 7. Fan (housing) | |

IMPORTANT NOTIFICATION

The 3 trays of the appliance shall be placed by respecting the indicated positions as on Picture.

A ferrite ring serving for EMC compatibility is mounted on the wire of the digital thermometer's sensor.

Do not remove the ferrite ring!

IN CASE IF YOUR REFRIGERATOR IS EQUIPPED WITH LOCK, DOMETIC Zrt DO NOT ACCEPT LIABILITY FOR LOSS OR DAMAGE OF KEYS. HOWEVER A COMPLETE LOCK CAN BE SUPPLIED.

3. Cleaning

i Clean the refrigerator thoroughly at intervals, as necessary. Remove the shelves; these, the cabinet interior and the door should be washed with a warm solution of sodium bi-carbonate. Finally wipe over with a cloth rinsed in warm water only, and dry with a clean cloth. Do not wash any plastic parts in water that is more than hand warm, and do not expose them to dry heat. The outside of the cabinet should be wiped down regularly with a clean, damp cloth.

NEVER USE STRONG CHEMICALS OR ABRASIVE CLEANING MATERIALS ON ANY PART OF THE CABINET.

4. Positioning the appliance

When first using this refrigerator please consider the following points:

- ENSURE REFRIGERATOR IS LEVEL
- THE PIPEWORK AT THE BACK GETS HOT IN USE
- THE REFRIGERATOR HAS NO 'MOTOR' AND IS THEREFORE VIRTUALLY SILENT
- ALLOW AT LEAST 3 HRS. FOR SIGNS OF COOLING

i During the refrigeration process, the appliance gives off heat from the condenser (under the upper part of the rear cover) into the surrounding air. The more ventilated the condenser is, the more effective the refrigeration will be.

- The other condition for satisfactory operation is that the appliance stands on a flat surface. This is best seen by placing a glass of water on top of the appliance.
- It is important that the appliance is not directly exposed to radiated heat (sunlight, radiator, near an oven, etc.).
- Please do not place the refrigerator onto thick carpet!

4.1. Installation instructions

This version is supposed to be freestanding. Cooling unit room cover protects from the hot parts of the cooling unit. Foot frame with grilles ensures sufficient ventilation.

5. Using the appliance

To start the refrigerator connect the plug to the socket outlet and switch on the socket. The cooling fins should show signs of cooling after about 1-2 hours depending on ambient temperature.

5.1. Initial operation

- The device must be plugged into a properly earthed socket.
- The miniCool starts automatically with a self-test. The LED illuminates after plugging-in several times and after that continuous light can be seen if the door is open.
- The miniCool achieves its operating temperature after several hours.

5.2. Temperature regulation

The Health Care refrigerator is fitted with a fully automatic electronic temperature regulator. It is not needed to set the temperature of the storage compartment manually.

6. Thermometer (Fig. 2)

Refer to the separate instruction leaflet supplied with the thermometer. The thermometer can be stood up using the attached stand or alternatively it can be fixed to the refrigerator door, in which case the sensor wire should be arranged so that it is not stretched or damaged when the door is opened.

The temperature sensor is located inside the refrigerator below the bottom shelf.

The display will then show the current, maximum and minimum temperature inside the refrigerator.

After the first pull down of the refrigerator (allow four hours) please make a "reset" on the thermometer because otherwise it will keep as "maximum" the ambient temperature.

The more medicines you store in the refrigerator then reduces the hazard of freezing, if the unit has a short period of time undercooling.

We take no responsibility for damage to vaccines through freezing, if there has been a cutout in the electricity for a short period of time.

The temperature readings may rise temporarily after opening the door or placing additional items inside the cabinet.

It is important to reset the thermometer regularly to ensure that the temperature is maintained between 2 °C and 8 °C.

On the battery transit strip must be removed to activate the temperature display.

Thermometer:

1. Min. Temperature
2. Max. Temperature
3. Current Temperature
4. Reset switch
5. Ferrite ring

7. Useful suggestion

The appliance is equipped with a fan in order to have a uniform inner temperature. The fan is operating only if the door is closed. At a door opening the fan will stop, then it starts again working when the door is closed.

If the fan incidentally stops, and the door of the appliance is closed, the electronics will emit an audible warning signal (continuous signal)

Audible warning signals to tell the customer to close the door: If the door is kept open for more than 20 s, the electronics emit an audible warning signal (intermittent signal with beeping frequency of 1 Hz), then after another 40 s it emits another audible warning signal (intermittent signal with beeping frequency of 2 Hz).

Do not overfill the appliance. It is advisable to leave some space between the products in the cabinet to ensure efficient and even cooling. If the shelves are packed too tightly there may be unacceptable temperature variations.

8. Defrosting, cleaning and maintenance (Fig. 3)

Frost will gradually form on the cooling surfaces and it must not be allowed to become too thick since it acts as an insulator and interferes with the cooling.

We recommend to defrost the appliance in every 3 weeks, or depending on the ice quantity on the cooling fins. While defrosting the appliance should be switched off by disconnecting the refrigerator from the electricity supply and leave the cabinet door open.

The frost will melt and the defrost water will run through the drain tube from the drip channel into an evaporating vessel located at the bottom rear of the refrigerator. (Fig 3.)

After defrosting the cabinet interior should be wiped out with a clean, damp cloth.

9. Interior light (Fig. 4)

Two light-emitting diodes (A) on an electronic PCB illuminate the cooling compartment when the door is opened.

Two sensors (B) control whether the door is open or not and the LEDs will work accordingly.

10. Replacing the decorative door panel (Fig. 5.)

If the refrigerator is furnished with handle, it is the handle that should be first removed at the change of the door panel.

Remove the mounter screws (1) of the forehead panel.(3)

Remove rosette (2) (in case of with lock version fridge). Lift and remove the forehead panel.(3)

Slide the decorative panel (4) out of the door frame and insert the new panel (thickness max. 2 mm).

Place back the forehead panel,rosette and screw in reverse sequence.

11. Fitting up and removing the handle (optional)

The handle which can be found among the accessories of the appliance should be fitted and removed as follows:

11.1. Fitting up the handle (fig. 6., 7.)

11.2. Removing the handle (fig. 8.)

12. Customer service

Before you call a service engineer, carry out the following simple tests:

- Is there a power failure? Check the fuses.
- Is the plug firmly connected to the socket and is the socket switched on? Check the socket by plugging in another appliance, e.g. a table lamp.
- Has the door not been closed properly causing the inside temperature to rise too quickly?
- Is the refrigerator level?
- Has the ventilation become obstructed?

If after checking all the above points, the refrigerator is still not working satisfactorily, contact your nearest service engineer. State the problem, model, product- and serial-number.

(Details of the model, product- and serialnumber are to be found on the data label in the top left of the inside compartment.)

This refrigerator is equipped with overheating protection. An authorised service technician can check whether this has been triggered or not.

In the event of a fault developing, contact your nearest Service Centre-see local telephone directory.

www.dometic.com

13. Putting into operation

Technical data:

Model	DS301H	DS601H
Type	DSH20-60	DSH20-60
Gross volume	28,1 litre	52,6 litre
Mains operation	220 - 240V (AC)	220 - 240V (AC)
Input	65W	85W
Energy consumption	0.86 kWh/24h	1.05 kWh/24h
Climate class	SN	SN
Refrigerant	151g H ₂ O + 101g NH ₃	157g H ₂ O + 105g NH ₃

13.1. Connecting to electrical mains

The appliance may only be operated from nominal voltage mains as shown on the data plate. The appliance mains plug may be connected to a mains socket earthed in accordance with regulations.

Any electrical work required to install this appliance should be carried out by a qualified electrician or competent person.

The manufacturer declines any liability should these safety measures not be observed.

Electrical Requirements

Before switching on, make sure the electricity supply voltage is the same as that indicated on the appliance rating plate.

Only for UK:

The appliance is supplied with a 3 amp plug fitted. In the event of having to change the fuse in the plug supplied, a 3 amp ASTA approved (BS 1362) fuse must be used. Should the plug need to be replaced for any reason, the wires in the mains lead are coloured in accordance with the following code:

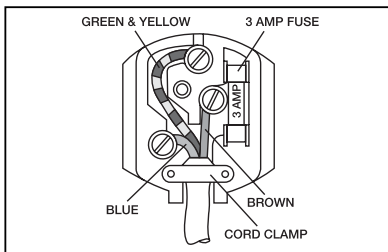
Green and Yellow:	Earth
Blue:	Neutral
Brown:	Live

The wire coloured green and yellow must be connected to the terminal marked with the letter „E“ or by the earth symbol \oplus or coloured green and yellow.

The wire coloured blue must be connected to the terminal „N“ or coloured black.

The wire coloured brown must be connected to the terminal marked „L“ or coloured red.

Upon completion there must be no cut, or stray strands of wire present and the cord clamp must be secure over the outer sheath.

14. Environmental protection information

The appliance does not contain any CFCs/HCFCs.

Ammonia (natural hydrogen and nitrogen compound) is used as a refrigerating agent in the refrigeration unit.

The ozone-friendly cyclopentane is activated as a motive agent for the PU foam insulation.

Sodium chromate is used for corrosion protection (less than 2 weight % of the coolant).

15. Recycling

After unpacking the appliance, the packing materials should be delivered to a local collection site. At the end of its useful lifetime, the appliance should be delivered to a specialised collection and reprocessing firm, which recclaims the usable materials. The rest is properly destroyed.

Appliances bearing this symbol must be deposited at the designated local reception point for the disposal of electrical and electronic equipment.

It is not permitted that this product be disposed of by way of the normal household refuse collection system.

Dometic refrigerators bear this symbol on the specifications plate (data plate) to be found in the interior of the unit.

This appliance complies with the following EEC directives:

LVD-Directive	2006/95/EC
EMC-Directive	2004/108/EC
CE-Directive	93/68/EEC
RoHS-Directive	2002/95/EC
WEEE-Directive	2002/96/EC

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